
# Food DNA Isolation Kit

**NORGEN**  
BIOTEK CORP.

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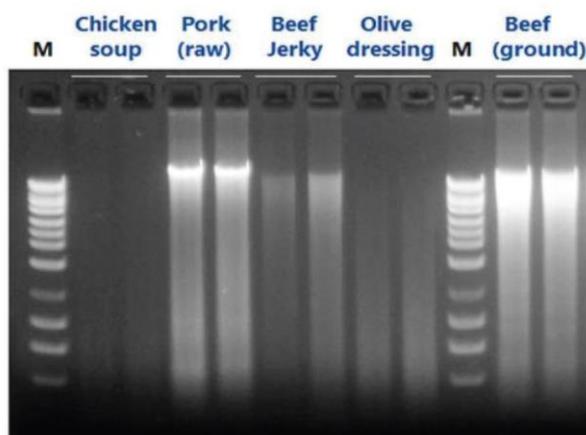
**Figure 1. Isolate DNA from a Wide Range of Food Samples.** Total genomic DNA was isolated from 200 mg of processed (chicken soup, beef jerky) or non-processed (raw pork, olive dressing, ground beef) food materials using Norgen's Food DNA Isolation kit. Following isolation, 10  $\mu$ L from each 100  $\mu$ L elution was loaded on 1% TAE agarose gel. Lane M: Norgen's HighRanger 1kb DNA Ladder.

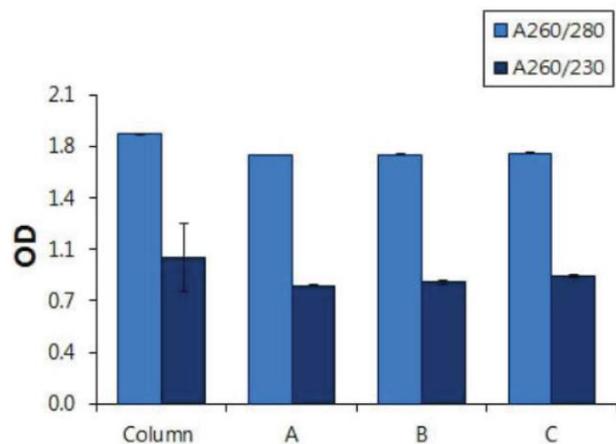**Figure 2. High Purity of DNA Samples.** DNA was isolated from 200 mg of processed (chicken soup, beef jerky) or non-processed (raw pork, olive dressing, ground beef) food materials using Norgen's Food DNA Isolation Kit. DNA purity was then determined using aNanoDrop. All the DNA isolated from the different food samples showed a high quality ( $260/280=1.9-2.1$ ).

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# Food DNA Isolation Kit

NORGEN  
BIOTEK CORP.

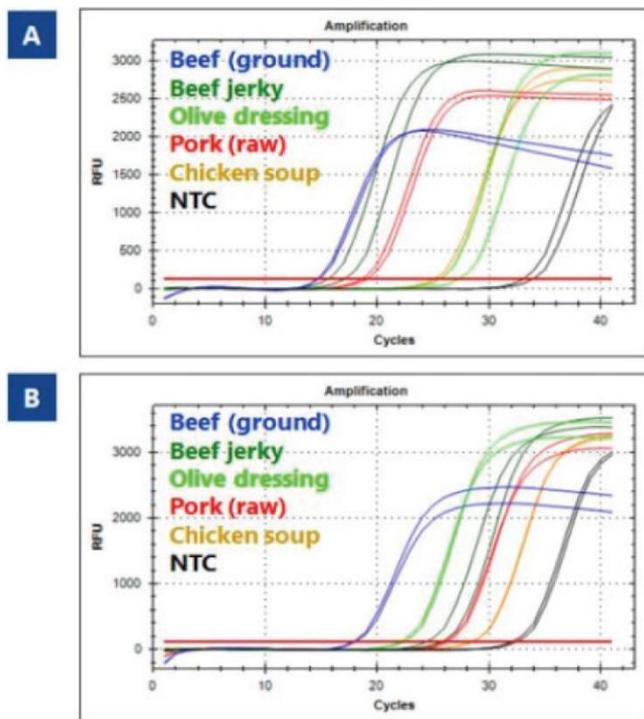
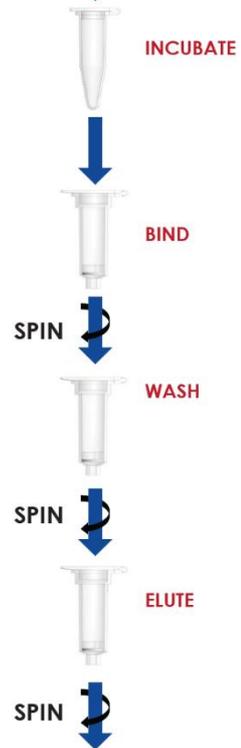
Figure 3. High Quality DNA free from PCR Inhibitors - Superior Quality to detect two house keeping genes (5S rDNA and 16S rDNA) representing GMO gene and food related pathogen detection. Total DNA was isolated from 200 mg samples of different food materials (processed and non-processed) using Norgen's Food DNA Isolation Kit. Next, 4  $\mu$ L of the elution was then used as the template in 20  $\mu$ L PCR reactions using universal 5s rDNA primers (A) and prokaryotic 16s rDNA primers (B) in real-time PCR (SYBR Green). Norgen's DNA was successfully amplified, indicating the high quality of the inhibitor-free DNA for GMO and pathogen detection from diversified food materials.

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